

Helpful Hints

1. Go to File. Select “IMPORT IMAGE”>”DIRECTORHOW CAPTURE” a video window is opened. You can also click on the video icon located at the top of the window.
2. Go to “Device” menu on the start page and click to select the appropriate model of the camera.
3. To capture video images:
 - a. Go to ‘CAPTURER” menu and select one of the following capture options,
 - i. Copy a frame (does not allow captured image to be saved)
 - ii. Capture a Frame (allows saving)
 - iii. Auto Capture
 - iv. Start Capture
 - b. To Capture a still image
 - i. Go to ‘FILE” select “IMPORT IMAGE”>’TWIN ACQUIRE”
4. You may need to go to “DEVICE” to select the correct eyepiece camera model.
5. You can select the file location and file type when saving by going to the “FILE” menu.