

Quick Start Cahn Microbalance

(revised 2016)

1. Turn on the balance with switch on rear
2. Let warm up 1 hour.
3. Choose the stirrup position to use
 - a. Stirrup position A on left end has max capacity of 250 mg
 - b. Stirrup position B close to center point has max capacity of 1250 mg
4. Choose sensitivity range or 1 and 0.1 μg for A or 10 μg for B.
5. Press the TARE button.
6. Place your sample on the stirrup.
7. When the weight display has stabilized read the weight.